



5X MORE
PRECISE
than standard
regulators



Research Series

$\frac{1}{16}$ " – $\frac{1}{4}$ " Back Pressure Regulators

LOW FLOW, LABORATORY SCALE FOR GAS, LIQUID, & MIXED PHASE SERVICE



The Equilibar Difference

OUR PERFORMANCE.

Equilibar® back pressure regulators outperform the competition particularly in applications with low flow rates, mixed phase fluids, corrosive media, or extreme temperatures.

OUR PEOPLE.

Every inquiry gets special attention from our Engineering team to determine the best possible product for your needs. Every back pressure regulator is hand assembled and tested to meet our stringent quality standards.

OUR PRIORITIES.


Our goal is to exceed your expectations. In an industry where delivery times frequently exceed 6 weeks, we offer many of our standard products with delivery in about a week.

Traditional back pressure regulators set the upstream pressure with a spring. These designs utilize sliding seals and other moving parts that can introduce hysteresis and other undesired effects into a process. The Equilibar® back pressure regulator uses a thin, supple diaphragm as the only moving part. This allows frictionless operation without cracking pressure or hysteresis. The accuracy of the Equilibar® back pressure regulator is determined by the accuracy of the pilot setpoint.




How it Works

Simply “load” the Equilibar® back pressure regulator with a pilot pressure equal to your desired back pressure and the Equilibar does the rest. This pressure forces the flexible diaphragm down onto a plate of orifices. A rise in inlet pressure lifts the diaphragm up to allow excess pressure to be relieved through the outlet orifices. Similarly, a loss of pressure at the inlet causes the diaphragm to be pushed closer to the orifices, restricting flow and rebuilding pressure upstream.

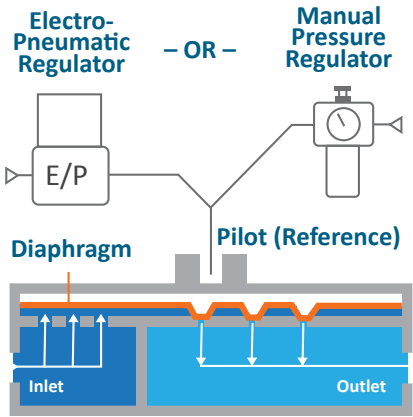

Pilot operate your Equilibar® back pressure regulator with an electronic pressure regulator for automated back pressure control.

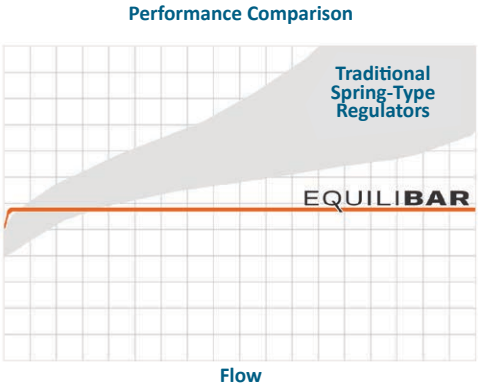

Or set the pilot pressure with a precision pressure reducing regulator for manual control.

| TYPE | PRESSURE REDUCING REGULATOR | BACK PRESSURE REGULATOR |
|-------------------|------------------------------|----------------------------|
| SCHEMATIC | | |
| CONTROLS PRESSURE | Downstream | Upstream |
| OPENS TO | Increase downstream pressure | Decrease upstream pressure |
| CLOSES TO | Decrease downstream pressure | Increase upstream pressure |

BACK PRESSURE REGULATORS VS PRESSURE REDUCING REGULATORS

Pressure reducing regulators reduce a higher supply pressure at the inlet down to a regulated lower pressure at the outlet (downstream). Back pressure regulators work the opposite way. They regulate the inlet (upstream) pressure by opening up only as much as necessary to hold back the desired pressure at the inlet (upstream).


Applications

There are hundreds of potential applications for the unique capabilities of Equilibar back pressure regulators. The Research Series is specifically designed for gas, liquid, and mixed phase applications where precision and consistency are critical.

Equilibar® Research Series regulators are especially useful in processes with low flow rates, extreme high pressures, and other challenging laboratory scenarios. By using unique combinations of diaphragm and O-ring materials, Equilibar regulators are able to perform in the harshest environments, including those with high temperatures and aggressive chemicals.


CATALYST RESEARCH REACTORS

Many catalyst research applications rely on high performance back pressure regulators to maintain stable pressure at elevated temperatures and down to nano flow rates. Because of the Equilibar's unique ability to handle two-phase flow, it is possible to substitute a low-pressure liquid/gas separator for the high-pressure separator.

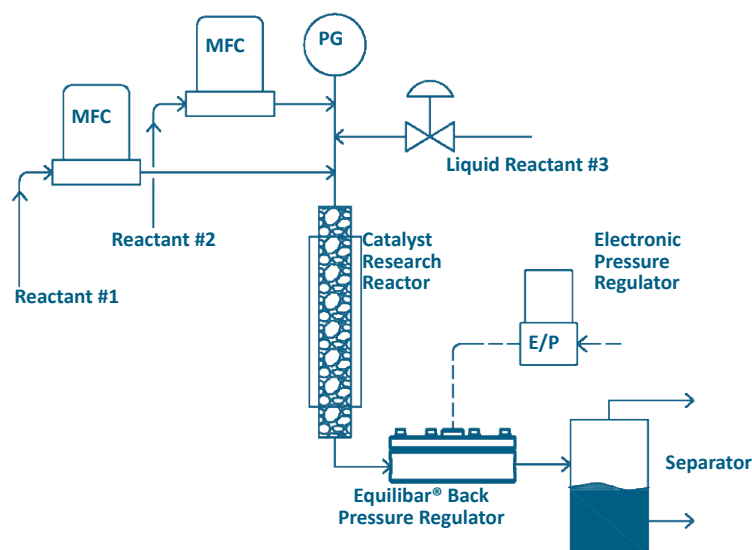

RESERVOIR CORE ANALYSIS

There are many types of upstream petroleum core analysis applications that depend on back pressure regulators to maintain the pressure of the sample. Core flooding (shown below), PVT analysis, pyrolysis, and slim tube analysis can all benefit from the ultra-low flow capability of Equilibar's Research Series. The ZF (Zero Flow) Series in particular was developed to address the demanding requirements of this industry.

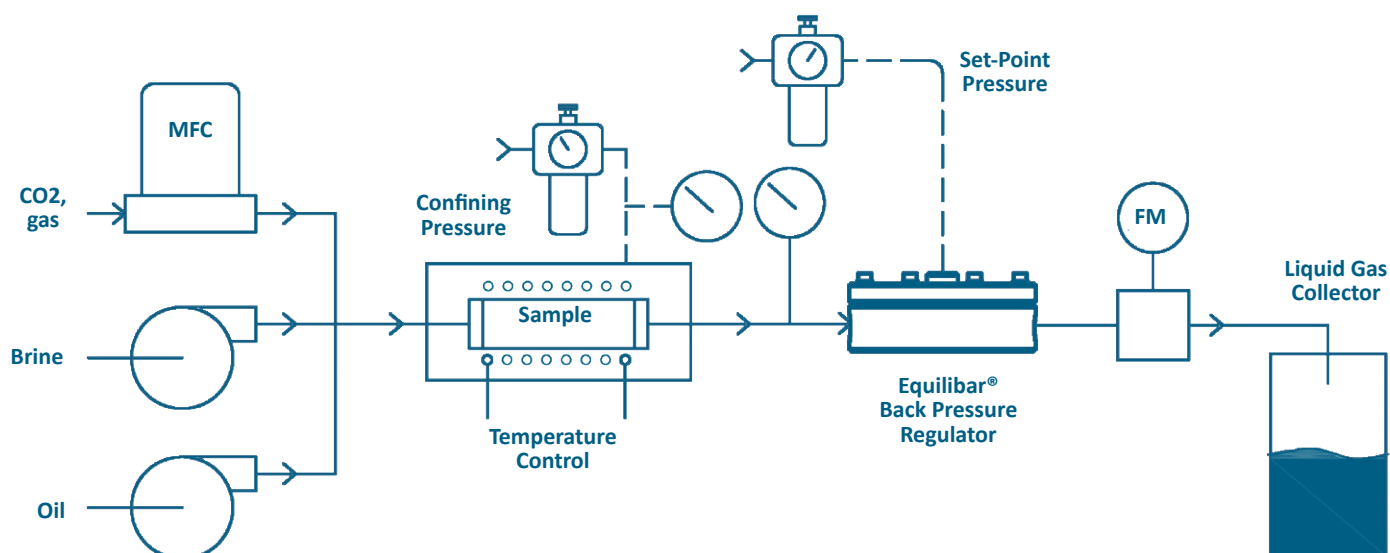

Key Advantages of Equilibar® Back Pressure Regulators

EXTREMELY WIDE RANGE OF CV

Control flow over 100,000:1 turndown ratio. Use just one Equilibar back pressure regulator to replace multiple control valves

HIGH PRESSURE CAPABILITIES

Equilibar Research Series back pressure regulators can operate at pressures up to 10,000 psig / 690 bar(g).

CHEMICAL COMPATIBILITY

Equilibar back pressure regulators are made in many exotic metal alloys and polymers including Hastelloy, Titanium, Zirconium, PTFE, PVDF, with dozens of diaphragm and O-ring options available.

HIGH TEMPERATURE CAPABILITIES

Equilibar Research Series back pressure regulators can be used up to 450C (840F) to allow hot media and/or oven installation.

MULTI-PHASE FLOW COMPATIBILITY

Traditional back pressure regulators consist of a single orifice and valve seat, while Equilibar back pressure regulators have multiple orifices in parallel. This patented multi-orifice design controls liquid and gas flow simultaneously without the pressure spikes seen in traditional regulators. This multi-orifice design allows pressure control of slurries and viscous fluids as well.

INSTANT RESPONSE TIME

Equilibar back pressure regulators instantly respond to changes in either line pressure or set point pressure.

Equilibar Research Series regulators are widely used in catalyst research systems because of their high temperature tolerance, two-phase flow capability, and ease of computer automation. (Photo courtesy Integrated Lab Solutions, GmbH)

RUGGED CONSTRUCTION & EASE OF SERVICE

Equilibar back pressure regulators are machined directly from bar stock metals.

Each Equilibar back pressure regulator consists of just a body, bolts, O-rings, and diaphragm.

Equilibar back pressure regulators can be completely cleaned and rebuilt in minutes. Rebuild kits are always in stock and ready to ship.

FRICTIONLESS OPERATION

Equilibar back pressure regulators are free of springs and valve seats that add friction and reduce precision.

LOW PRESSURE CAPABILITIES


Equilibar's direct diaphragm sealing technology controls down to 0.5 in WC.

COMPUTER AUTOMATION

All Equilibar back pressure regulators can be controlled both manually and electronically. Verify the system design with manual control, then automate it later by adding an electronic pilot regulator

DIRECT CUSTOMER SUPPORT

Contact an [Equilibar application engineer](#) to assist with questions or issues.

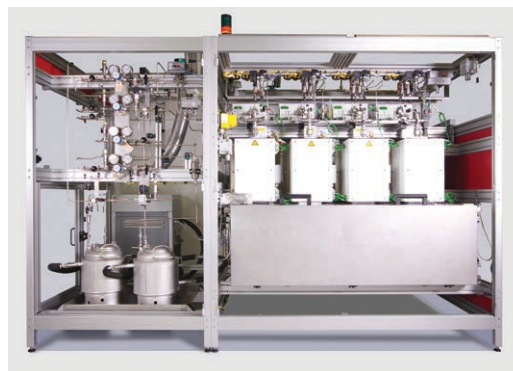

Standard Multi-Orifice Series Back Pressure Regulators

Our unique, patented multi-orifice technology has an extremely wide flow range and superior two phase flow performance.

| BASE | MAX PRESSURE RATING | FLOW COEFF. (CV) | | INLET/OUTLET PORT SIZE | REFERENCE PORT SIZE | PORT THREADS | | DIM A | DIM B | WEIGHT | Dead Volume ¹ |
|--------------------------|---------------------|------------------|------|---------------------------|------------------------|--------------|---------------------------|-----------|----------|-----------|--------------------------|
| PART # | PSIG (BAR) | MIN | MAX | | | STANDARD | | INCH (MM) | | LB (KG) | mL (in ³) |
| STANDARD LOW FLOW MODELS | | | | | | | | | | | |
| LF0 | 1000 (68) | 1E-08 | 0.01 | 1/16'' | 1/8" | V (HPLC) | A, C, N | 2.5 (64) | 1.5 (39) | 1.7 (0.8) | 0.23 (0.014) |
| LF1 | | 1E-08 | 0.07 | 1/8" | | N (NPT) | A, B, C, O, R, T, V, W | 2.5 (64) | 1.5 (39) | 1.7 (0.8) | 2.19 (0.134) |
| LF2 | | 1E-08 | 0.07 | 1/4" | | | | 2.5 (64) | 1.5 (39) | 1.7 (0.8) | 3.27 (0.200) |
| H3P0 | 3000 (200) | 1E-08 | 0.01 | 1/16'' | 1/8" | V (HPLC) | A, C, N | 2.5 (64) | 1.7 (42) | 2.0 (0.9) | 0.17 (0.010) |
| H3P1 | | 1E-08 | 0.07 | 1/8" | | N (NPT) | A, B, C, O, R, T, V, W | 2.5 (64) | 1.7 (42) | 2.0 (0.9) | 2.08 (0.127) |
| H3P2 | | 1E-08 | 0.07 | 1/4" | | | | 2.8 (70) | 1.7 (42) | 2.5 (1.1) | 3.48 (0.212) |
| H6P0 | 6000 (400) | 1E-08 | 0.01 | 1/16'' | 1/8" | V (HPLC) | A, C, N | 2.8 (70) | 1.7 (42) | 2.5 (1.1) | 0.17 (0.010) |
| H6P1 | | 1E-08 | 0.07 | 1/8" | | N (NPT) | A, B, C, O, R, T, V, W | 2.8 (70) | 1.7 (42) | 2.5 (1.1) | 2.18 (0.133) |
| H6P2 | | 1E-08 | 0.07 | 1/4" | | | | 2.8 (70) | 1.7 (42) | 2.5 (1.1) | 3.25 (0.198) |
| H10P1 | 10,000 (680) | 1E-06 | 0.07 | 1/8" | 1/8" | W | A, C | 3.0 (76) | 2.1 (53) | 3.0 (1.4) | 1.28 (0.078) |
| HIGHER FLOW MODELS | | | | | | | | | | | |
| HF1 | 1000 (68) | 1E-05 | 0.41 | 1/8" | 1/8" | N (NPT) | A, B, C, O, R, T, V, W | 2.5 (64) | 1.5 (39) | 1.7 (0.8) | 5.65 (0.345) |
| HF2 | | | 0.41 | 1/4" | | | | 2.5 (64) | 1.5 (39) | 1.7 (0.8) | 6.68 (0.408) |
| H3PF2 | 3000 (200) | 1E-05 | 0.35 | 1/4" | 1/8" | N (NPT) | A, B, C, O, R, T, V, W | 2.8 (70) | 1.7 (42) | 2.5 (1.1) | 7.75 (0.473) |
| H6PF2 | 6000 (400) | | 0.35 | 1/4" | | | | 2.8 (70) | 1.7 (42) | 2.5 (1.1) | 6.23 (0.380) |
| HIGH TEMPERATURE MODELS | | | | | | | | | | | |
| HT1 | 5000 (345) | 1E-05 | 0.07 | 1/8" | 1/8" | N (NPT) | A, B, C, O, R, T, V, W | 3.3 (82) | 2.0 (50) | 3.5 (1.6) | 2.62 (0.160) |
| HT2 | | 1E-05 | 0.07 | 1/4" | | | | 3.8 (95) | 2.0 (50) | 4.3 (2.0) | 3.70 (0.226) |
| HTF1 | 5000 (345) | 1E-05 | 0.35 | 1/8" | 1/8" | N (NPT) | A, B, C, O, R, T, V, W | 3.8 (95) | 2.0 (50) | 4.3 (2.0) | 8.93 (0.545) |
| HTF2 | | 1E-05 | 0.35 | 1/4" | | | | 3.8 (95) | 2.0 (50) | 4.3 (2.0) | 8.93 (0.545) |

¹ Dead volumes are approximate and for reference only. Approximations are made assuming diaphragm is in the "down" position.

| PORTING OPTIONS | | |
|-----------------|----------------------|--------------|
| Notation | Type | Max Cv |
| N | NPT (Standard) | Full |
| A | HiP (High Pressure) | 0.072 |
| B | BSPP | Full |
| C | Custom | — |
| O | Swagelok VCO* | 0.072 |
| R | Swagelok VCR* | 0.072 |
| T | Tube Stub | 0.072 |
| V | HPLC | .001 - 0.018 |
| W | Autoclave Speed-Bite | .07 |

| TECHNICAL SPECIFICATIONS | |
|--------------------------|---|
| Max Operating Pressure | Pressure ratings listed in the table are the maximum possible pressure that a unit may be configured to. Units can be configured for optimum performance at lower pressures. Speak with an Application Engineer for more information. |
| Proof Pressure | 150% Rated Pressure ¹ |
| Design Pressure | 400% Maximum Body Pressure ² |
| Temperature Capabilities | Up to 150C (Metallic Body, PTFE Diaphragm, Viton® O-rings) Up to 200C (Metallic Body, Metallic Diaphragm, Viton® O-rings) Up to 300C (Metallic Body, Metallic Diaphragm, Kalrez® O-rings) Up to 450C (HT/HTF Models, Metallic Body, Grafoil Gaskets) |


¹All Equilibr units are tested to 150% of their rated pressure prior to shipment.

²Designed according to ASME B31.3, which incorporates an approximate 4X safety factor.

³Polymer units reduce maximum allowable working pressure; consult an application engineer for additional information.

Viton® and Kalrez® are registered trademarks of DuPont. VCO® and VCR® are trademarks of Swagelok. Grafoil® is a trademark of GrafTech.


⚠ Equilibr regulators are control devices, not safety devices or shut-off devices and should not be used as such.

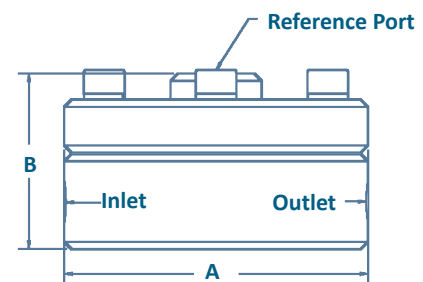


Mounting brackets available for most models. Consult an Application Engineer for mounting information.

| WETTED MATERIALS | |
|------------------|--|
| Body Material | Stainless Steel 316/316L (standard) Also available: Hastelloy C276, Titanium, Zirconium, PTFE ³ , PVDF ³ , PEEK ³ , PVC ³ , Monel |
| O-Rings | Viton® (FKM) (standard) Also available: Kalrez® (FFKM), PTFE, EPDM, Buna-N, Grafoil® (HT Only) |
| Diaphragm | PTFE/Glass Laminate (standard) Also available: Stainless Steel SS316/316L, Hastelloy C276, Virgin PTFE, FKM, Polyimide, Buna-N, PEEK, EPDM |

PERFORMANCE¹ OF STANDARD LOW FLOW MODELS (NITROGEN)

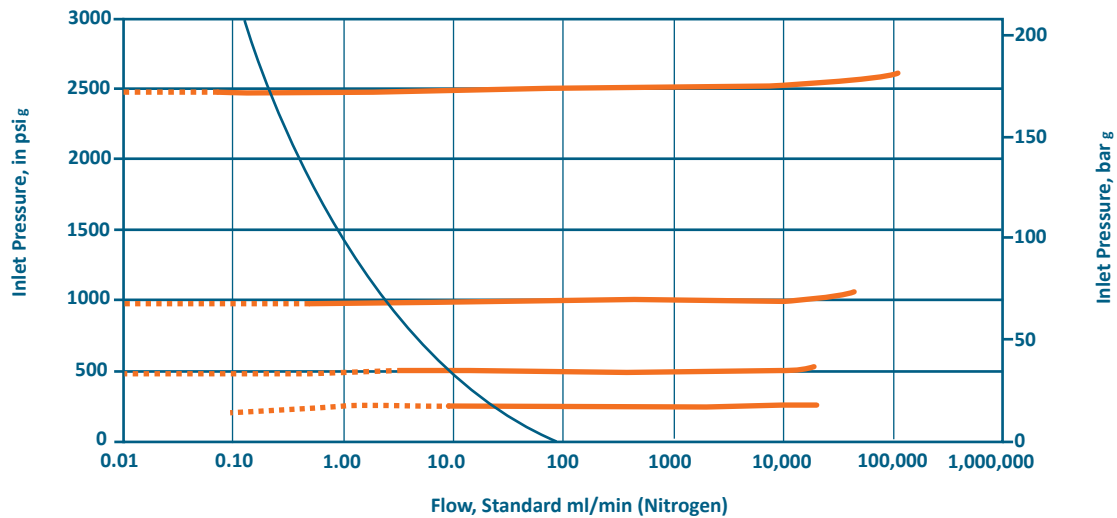

¹Performance below and left of the blue curve is configuration dependent. Higher flow models exhibit similar performance at greater flow capacity. While these performance curves are for nitrogen, Equilibar Back Pressure Regulators may be used for gas, liquid, or mixed phase service.

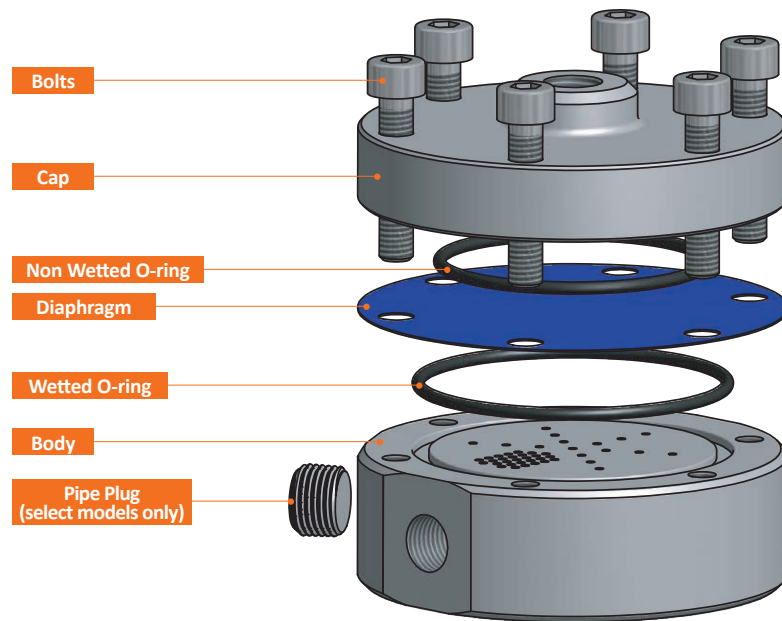

Ultra Low Flow Series Back Pressure Regulators

The Ultra Low Flow Series performs similarly to the Standard Multi-Orifice Series but can operate at lower flow rates.

| BASE | MAX PRESSURE RATING | FLOW COEFF. (CV) | | INLET/OUTLET PORT SIZE | REFERENCE PORT SIZE | PORT THREADS | | DIM A | DIM B | WEIGHT | Dead Volume ¹ |
|--------------------------------------|---------------------|------------------|------|---------------------------|------------------------|--------------|---------------------------|-----------|----------|-----------|--------------------------|
| PART # | PSIG (BAR) | MIN | MAX | | | STANDARD | OPTIONAL | INCH (MM) | | LB (KG) | mL (in³) |
| U3L0 | 3000 (200) | 1E-09 | 0.01 | 1/16" | 1/8" | V (HPLC) | A, N, C | 2.5 (64) | 1.7 (42) | 2.0 (0.9) | 0.25 (0.015) |
| U3L1 | | 1E-09 | 0.05 | 1/8" | | N (NPT) | A, B, C, O, R, T, V, W | 2.5 (64) | 1.7 (42) | 2.0 (0.9) | 2.16 (0.132) |
| U3L2 | | 1E-09 | 0.05 | 1/4" | | | | 2.8 (70) | 1.7 (42) | 2.5 (1.1) | 3.34 (0.204) |
| U6L0 | 6000 (400) | 1E-09 | 0.01 | 1/16" | 1/8" | V (HPLC) | A, N, C | 2.8 (70) | 1.7 (42) | 2.5 (1.1) | 0.25 (0.015) |
| U6L1 | | 1E-09 | 0.05 | 1/8" | | N (NPT) | A, B, C, O, R, T, V, W | 2.8 (70) | 1.7 (42) | 2.5 (1.1) | 2.28 (0.139) |
| U6L2 | | 1E-09 | 0.05 | 1/4" | | | | 2.8 (70) | 1.7 (42) | 2.5 (1.1) | 3.34 (0.204) |
| U10L0 | 10,000 (680) | 1E-06 | 0.01 | 1/16" | 1/8" | V (HPLC) | A, N, C | 3.0 (76) | 2.1 (53) | 3.0 (1.4) | 3.34 (0.204) |
| U10L1 | | 1E-06 | 0.05 | 1/8" | | N (NPT) | A, C, W | 3.0 (76) | 2.1 (53) | 3.0 (1.4) | 1.33 (0.081) |
| HIGH TEMPERATURE MODELS ² | | | | | | | | | | | |
| ULHT1 | 5000 (345) | 1E-09 | 0.05 | 1/8" | 1/8" | N (NPT) | A, B, C, O, R, T, V, W | 3.3 (84) | 2.0 (51) | 4.5 (2.0) | 2.44 (.149) |
| ULHT2 | | 1E-09 | 0.05 | 1/4" | | | | 3.8 (96) | 2.0 (51) | 5.2 (2.4) | 3.54 (.216) |

¹ Dead volumes are approximate and for reference only. Approximations are made assuming diaphragm is in the "down" position.

² ULHT models are for ultra low flow / high temperature applications up to 450C with specific gaskets and diaphragm.


PORTING OPTIONS

| Notation | Type | Max Cv |
|----------|---------------------------|--------------|
| N | NPT (Standard) | Full |
| A | HiP (High Pressure) | Full |
| B | BSPP | Full |
| C | Custom | — |
| O | Swagelok VCO* | Full |
| R | Swagelok VCR* | Full |
| T | Tube Stub | Full |
| V | HPLC | .001 - 0.018 |
| W | Autoclave Speed-Bite W125 | Full |

TECHNICAL SPECIFICATIONS


| | |
|------------------------|---|
| Max Operating Pressure | Pressure ratings listed in the table are the maximum possible pressure that a unit may be configured to. Units can be configured for optimum performance at lower pressures. Speak with an Application Engineer for more information. |
| Proof Pressure | 150% Rated Pressure ¹ |
| Design Pressure | 400% Maximum Body Pressure ² |
| Temperature Rating | 150C (Metallic Body, PTFE Diaphragm, Viton® O-rings) 200C (Metallic Body, Metallic Diaphragm, Viton® O-rings) 300C (Metallic Body, Metallic Diaphragm, Kalrez® O-rings) Up to 450C (ULHT Models, Metallic Body, Grafoil Gaskets) |

Performance⁴ of Ultra Low Flow Series (Nitrogen)

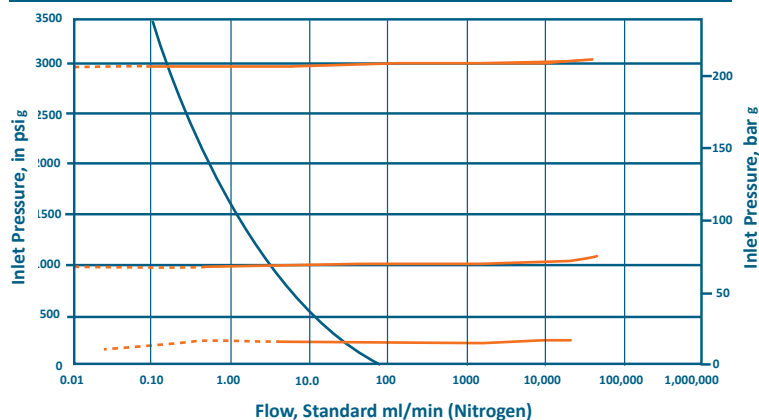

⁴Performance below and left of the blue curve is configuration dependent. While the above curves are for nitrogen, Equilibrar Back Pressure Regulators may be used for gas, liquid, or mixed phase fluids.

Reference Port

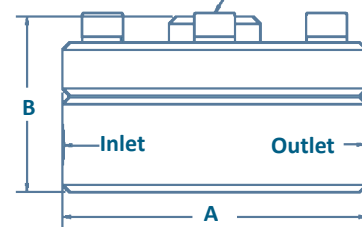

Mounting brackets available for most models. Consult an application engineer for mounting information.

WETTED MATERIALS

| | |
|---------------|--|
| Body Material | Stainless Steel 316/316L (standard) Also available: Hastelloy C276, Titanium, Zirconium, PTFE ³ , PVDF ³ , PEEK ³ , PVC ³ , Monel |
| O-Rings | Viton® (FKM) (standard) Also available: Kalrez® (FFKM), PTFE, EPDM, Buna-N |
| Diaphragm | PTFE/Glass Laminate (standard) Also available: Stainless Steel SS316/316L, Hastelloy C276, Virgin PTFE, FKM, Polyimide, Buna-N, PEEK, EPDM |


¹All Equilibrar units are tested to 150% of their rated pressure prior to shipment.

²Designed according to ASME B31.3, which incorporates an approximate 4X safety factor.

³Polymer units reduce maximum allowable working pressure, consult an application engineer for additional information.

Viton® and Kalrez® are registered trademarks of DuPont. VCO* and VCR* are trademarks of Swagelok.

⚠ Equilibrar regulators are control devices, not safety devices or shut-off devices and should not be used as such.

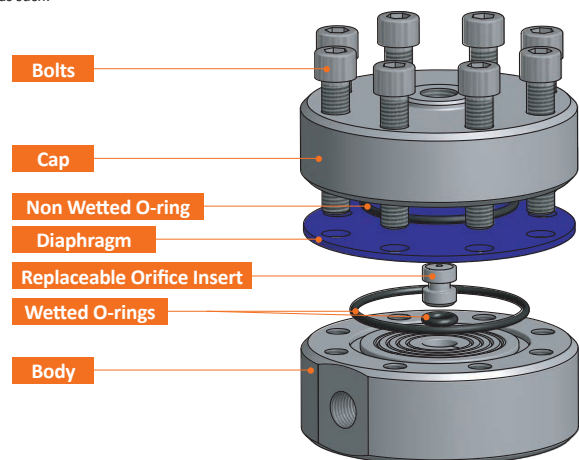

Zero Flow Series Back Pressure Regulator

Operates similarly to the Standard Multi-Orifice design, but utilizes a soft-seat to provide effective zero flow pressure control.

| BASE | MAX PRESSURE RATING | FLOW COEFF. (CV) | | INLET/OUTLET PORT SIZE | REFERENCE PORT SIZE | PORT THREADS | | DIM A | DIM B | WEIGHT | Dead Volume ¹ |
|--------|---------------------|------------------|-------|------------------------|---------------------|--------------|------------------------|-----------|----------|-----------|--------------------------|
| PART # | PSIG (BAR) | MIN | MAX | | | STANDARD | OPTIONAL | INCH (MM) | | LB (KG) | mL (in ³) |
| ZF0 | 3000 (200) | 0 | 0.015 | 1/16" | 1/8" | V (HPLC) | A, C, N | 2.5 (64) | 1.7 (42) | 2.0 (0.9) | 0.25 (0.02) |
| ZF1 | 3000 (200) | 0 | | 1/8" | 1/8" | N (NPT) | A, B, C, O, R, T, V, W | 2.5 (64) | 1.7 (42) | 2.0 (0.9) | 2.16 (0.13) |

¹ Dead volumes are approximate and for reference only. Approximations are made assuming diaphragm is in the "down" position. Speak with an Application Engineer to find out if the Zero Flow Series is right for your application.

| PORTING OPTIONS | | |
|-----------------|---------------------------|--------------|
| Notation | Type | Max Cv |
| N | NPT (Standard) | Full |
| A | HiP (High Pressure) | Full |
| B | BSPP | Full |
| C | Custom | — |
| O | Swagelok VCO [®] | Full |
| R | Swagelok VCR [®] | Full |
| T | Tube Stub | Full |
| V | HPLC | .001 - 0.018 |
| W | Autoclave Speed-Bite W125 | Full |

| TECHNICAL SPECIFICATIONS | |
|--------------------------|---|
| Max Operating Pressure | Pressure ratings listed in the table are the maximum possible pressure that a unit may be configured to. Units can be configured for optimum performance at lower pressures. Speak with an Application Engineer for more information. |
| Proof Pressure | 150% Rated Pressure ¹ |
| Design Pressure | 400% Maximum Body Pressure ² |
| Temperature Rating | 200C (Metallic Body, Metallic Diaphragm, Viton [®] O-rings) 300C (Metallic Body, Metallic Diaphragm, Kalrez [®] O-rings) |

| WETTED MATERIALS | |
|------------------|--|
| Body Material | Stainless Steel 316/316L (standard) Also available: Hastelloy C276, Titanium, Zirconium, PTFE ³ , PVDF ³ , PEEK ³ , PVC ³ , Monel |
| O-Rings | Viton [®] (FKM) (standard) Also available: Kalrez [®] (FFKM), PTFE, EPDM, Buna-N (Nitrile) |
| Diaphragm | Stainless Steel SS316/316L (standard) Also available: Hastelloy C276, PTFE/Glass Laminate, Virgin PTFE, FKM, Polyimide, Buna-N, (Nitrile) PEEK, EPDM |

¹All Equilibar units are tested to 150% of their rated pressure prior to shipment.


²Designed according to ASME B31.3, which incorporates an approximate 4X safety factor.

³Polymer units reduce maximum allowable working pressure, consult an application engineer for additional information.

Viton[®] and Kalrez[®] are registered trademarks of DuPont. VCO[®] and VCR[®] are trademarks of Swagelok.

⚠ Equilibar regulators are control devices, not safety devices or shut-off devices and should not be used as such.


Performance of Zero Flow Series (Nitrogen)⁴

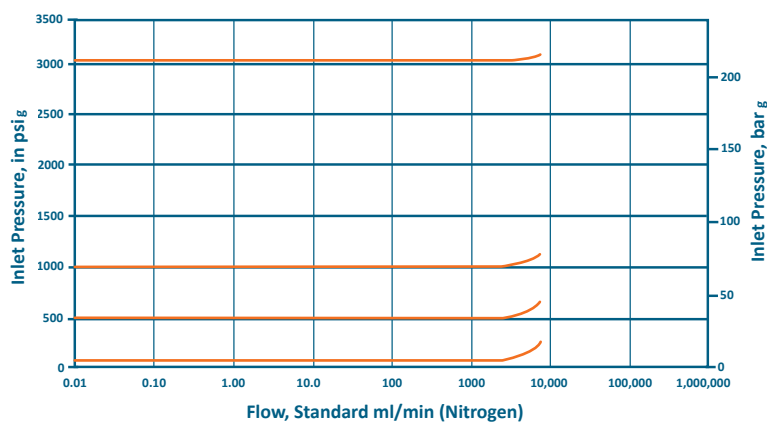


⁴While these performance curves are for nitrogen, Equilibar Back Pressure Regulators may be used for gas, liquid, or mixed phase service.

PATENTS

These regulators are subject to one or more of [these patents](#): US6,886,591, US7,080,660, US7,673,650, US8,215,336, US9,447,890 DE60322443D1, GB1639282, FR1639282, EP2724060.

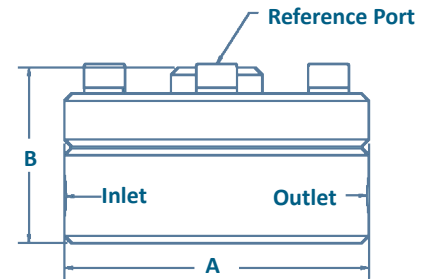

Mounting brackets available for most models. Consult an Application Engineer for mounting information.

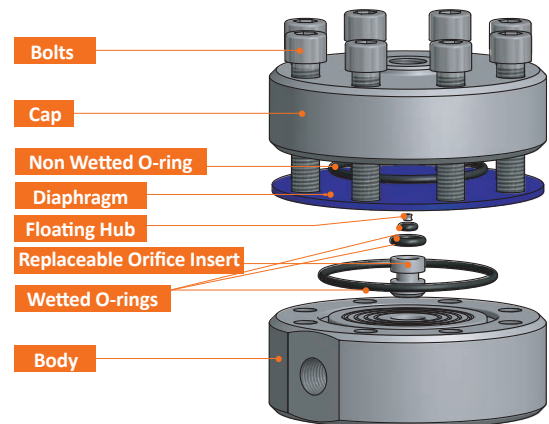

Research Series Part Number Key

This part number key explains our part numbering system and possible model options. All of our BPRs are custom-configured by our engineers based on the customer's specific application's parameters (process fluid, pressures, flow rates, temperature, etc.). Our engineers will request process operating parameters in order to build and quote a full part number for a suitable regulator. This chart is a reference to help understand the chosen part number

| EXAMPLE | | | | | | | | | | | | | | | | | | | |
|---------|---|---|---|---|---|---|---|---|---|---|-----|---|-----|----|----|----|----|----|----|
| LF | 1 | S | N | N | X | — | N | S | X | P | 500 | T | 150 | G | X | V | V | | B |
| | | | | N | X | — | | | X | P | | T | | | X | | | | |
| 1 | 2 | 3 | 4 | 5 | 6 | — | 7 | 8 | 9 | | 10 | | 11 | 12 | 13 | 14 | 15 | 16 | 17 |

| | | |
|--|---|---|
| <p>1 Model Up to 1000 psig LF Cv: 1E-8 to 0.07 HF Cv: 1E-5 to 0.41</p> <p>Up to 3000 psig U3L Cv: 1E-9 to 0.05 H3P Cv: 1E-8 to 0.07 H3PF Cv: 1E-5 to 0.35</p> <p>Up to 6000 psig U6L Cv: 1E-9 to 0.05 H6P Cv: 1E-8 to 0.07 H6PF Cv: 1E-5 to 0.35</p> <p>Up to 10,000 psig U10L Cv: 1E-6 to 0.05 H10P Cv: 1E-6 to 0.07 H10PF Cv: 1E-4 to 0.35</p> <p>Specialty Regulators HT High Temperature ZF Zero-Flow</p> <p>2 Port Size 0 1/16" 1 1/8" 2 1/4"</p> <p>3 Body Material S Stainless Steel 316/316L H Hastelloy C276 T Titanium Z Zirconium F PTFE K PEEK D PVDF</p> <p>4 Port Threads N NPT B BSPP W Autoclave Speed-Bite O VCO® R VCR® V HPLC A HiP</p> <p>5 Recess N None</p> | <p>6 Mod # X (Factory Selected)</p> <p>7 Reference Port Threads N NPT B BSPP W Autoclave Speed-Bite O VCO® R VCR® V HPLC A HiP</p> <p>8 Cap Material (Non Wetted) S Stainless Steel 316/316L P PVC F PTFE K PEEK D PVDF</p> <p>9 Bolts X (Factory Selected)</p> <p>10 Pressure Rating (psi_g) This is the maximum pressure you would like your unit to be configured to accept. Must be equal to or less than the maximum body pressure (in psi_g)</p> <p>11 Temperature Rating Temperature Limitations: 40C for most polymer bodies 150C for most PTFE diaphragms 200C for Viton O-rings 300C for Kalrez O-rings</p> <p>12 Diaphragm Material G PTFE (Glass Reinforced) B Buna-N (Nitrile) V FKM Fluoroelastomer M EPDM E Polyethylene F PTFE (Virgin) S Stainless Steel 316/316L H Hastelloy C276 I Polyimide K PEEK L Kel-F C FEP</p> <p>13 Diaphragm Thickness (Factory Selected)</p> | <p>14 O-ring (Wetted) V Viton® FKM Shore 75 W Viton® FKM Shore 90 K Kalrez® FFKM Grade 7075 L Kalrez® FFKM Grade 7090 F PTFE Grade 55 E EPDM Grade 70 B Buna-N Grade 70 (Nitrile) G Grafoil (HT Only)</p> <p>15 O-ring (Non Wetted) V Viton® FKM Shore 75 W Viton® FKM Shore 90 K Kalrez® FFKM Grade 7075 L Kalrez® FFKM Grade 7090 F PTFE Grade 55 E EPDM Grade 70 B Buna-N Grade 70 (Nitrile) G Grafoil® (HT Only)</p> <p>16 O-ring (Wetted) U3L, U6L, U10L, ZF Only Leave blank for all other models (Blank) None V Viton® FKM Shore 75 W Viton® FKM Shore 90 K Kalrez® FFKM Grade 7075 L Kalrez® FFKM Grade 7090 E EPDM Grade 70 B Buna-N Grade 70 (Nitrile)</p> <p>17 O-ring (Wetted) ZF Only Leave blank for all other models (Blank) None V Viton® FKM Shore 75 W Viton® FKM Shore 90 K Kalrez® FFKM Grade 7075 L Kalrez® FFKM Grade 7090 E EPDM Grade 70 B Buna-N Grade 70 (Nitrile)</p> <p>18 Special Options (Blank) None B Mounting Bracket O Oxygen Cleaning</p> |
|--|---|---|

Options listed in blue are typically in stock for quick shipment.

For custom options [contact an applications engineer](#).

Viton® and Kalrez® are registered trademarks of DuPont. VCO® and VCR® are trademarks of Swagelok. Grafoil® is a trademark of GrafTech.

About Equilibar

Equilibar provides innovative and robust pressure control technology for researchers and engineers worldwide. We are proud to design, manufacture and test our patented back pressure regulators in our factory overlooking the Blue Ridge Mountains near Asheville, NC.

APPLICATION ENGINEERING – WHAT SETS US APART

Unlike mass-market regulator distributors, we focus on working with you, the scientist or engineer with a complex pressure control scenario.


Our application engineers work collaboratively with clients to identify the optimal model, trim, and diaphragm for each application's unique challenges. No matter where you are on the globe, you can stay in close contact with your engineer by email, telephone, video-conference or fax.

After installation, your application engineer will support you with start-up information and fine-tuning as needed.

Equilibar, LLC
320 Rutledge Rd.
Fletcher, North Carolina 28732
United States
Tel: +1-828-650-6590
Fax: +1-801-504-4439
Monday - Friday
8:00 AM - 5:00 PM EST
12:00 - 21:00 GMT
inquiry@equilibar.com



Each application is reviewed by our engineering team to ensure quality performance of our products.

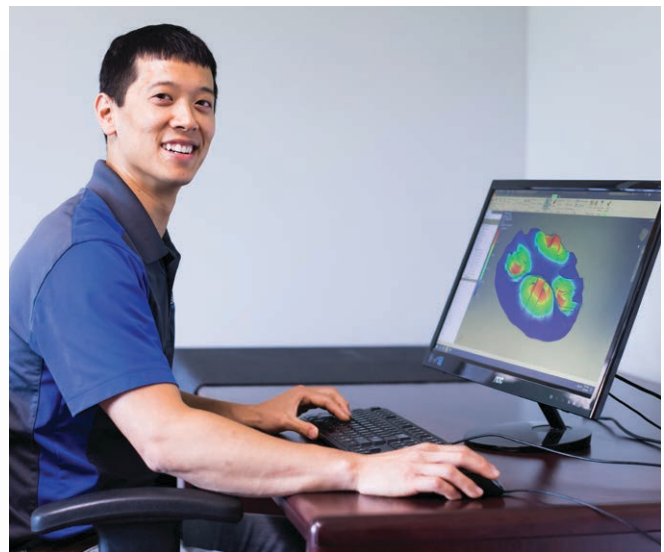

Our engineers offer custom designed solutions for the most difficult pressure control challenges. Feel free to contact us to discuss your situation.



320 Rutledge Rd. • Fletcher, NC 28732 • Tel: (828)650-6590 • Fax: (801)504-4439 • www.equilibar.com



Made in the
USA

Equilibar's quality system is
ISO 9001:2015 certified.